


## Using the 600ml beaker and beaker stand

The tank should be at least half filled with fluid. It is most important that the fluid is NOT water alone, but water and detergent or another cleaning fluid. The stand should be placed on top of the unit with the locating lugs pointing down. Beaker stands are available in single and double beaker formats. The rubber ring should be slid up or down the beaker so that when put into the beaker stand the base of the beaker is between 2 1/2cm (1") and 5cm (2") below the surface of the fluid but not touching the base of the tank. CARE SHOULD BE TAKEN not to lower the beaker too far as the fluid may flow over the top of the tank. The ultrasonic cleaner may then be operated.

This method is particularly useful for cleaning small components.

The single beaker stand and the basket can normally be used simultaneously\*

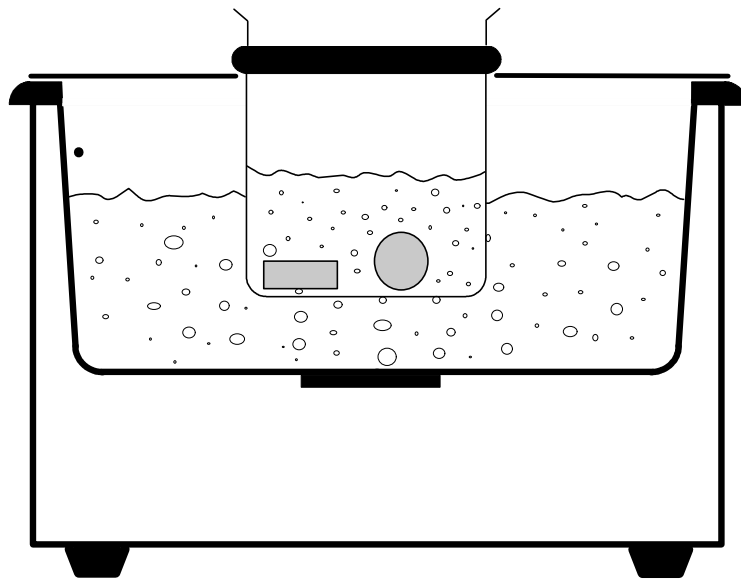

\*Dependant on basket design.